


appleton®

For the Life Scientist

Appleton appROCK 2D Rocker

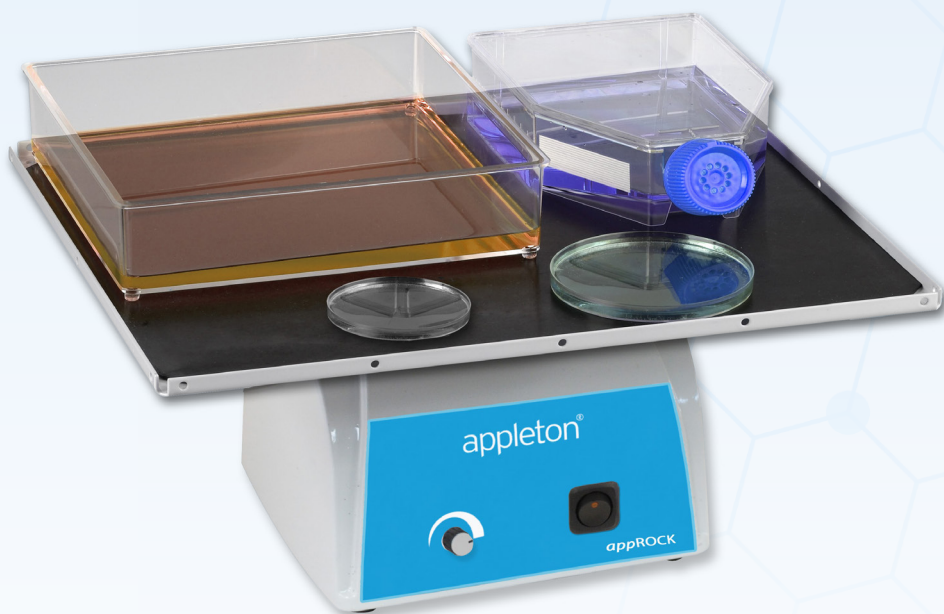

Instruction Manual

Product code: AER0001

Product Specifications

Speed:	Variable 2 to 30 rpm
Tilt Angle:	Variable 0 to +/-30°C
Operating Temperature:	+4 to 40°C
Platform:	35 x 30 cm
Load Capacity:	2.2 kg
Dimensions:	35 x 30 x 20cm
Weight:	3.6 kg
Electrical:	230V, 50 Hz
Warranty:	2 years

Product Set-Up

Place the rocker on a clean, flat, stable surface. Make sure that the power switch is in the OFF (not illuminated) position. Set the unit to the desired tilt angle (see below). Plug the rocker into a properly grounded outlet. Check that nothing blocks the up and down rocking action of the platform. **Set the Speed Control knob to the lowest speed setting (full counter clockwise).**

Product Operation

- Load the platform from the centre out, making sure the items placed on the platform create a balanced load around the centre line of the platform.
- Turn the unit on by activating the ON/OFF switch so that it is illuminated.
- Rotate the speed control knob clockwise to set the desired rocking speed.

Tilt Angle Adjustment

The tilt angle is adjusted at the back of the unit. First loosen the Safety Shield knob and move the Safety Shield aside to gain access to the tilt adjustment screw. Loosen the tilt adjustment screw with the 2.5 mm hex Allen wrench provided with the unit. Sliding the black adjustment bar will increase or decrease the platform tilt angle. Once the platform angle is selected, re-tighten the tilt angle adjustment screw with the Allen wrench. **VERY IMPORTANT!** Position the safety shield so that it is centred over the black oval adjustment bar, then hand tighten the Safety Shield knob. The safety shield is used to keep fingers out of the rotating mechanism.

Accessories, stacking platforms and mats

A stacking platform is available, which effectively doubles the work area. There is a 7.5 cm separation between the two platforms, allowing access to the content of both trays. The stacking platform comes with a flat mat. Please contact us for more information.

Care and Maintenance

No routine maintenance is required other than to keep the unit clean. Cleaning can be done with a damp cloth. Avoid the use of harsh solvents. Appleton Woods distributes these products in the UK, for customer service or technical support, please contact: info@appletonwoods.co.uk or call on **0121 4587740**.

Warranty

This product is covered by the **Appleton Promise** which means that should any faults arise within the 2 years warranty period it will be replaced free of charge. Damage from neglect, misuse or dropping the unit will not be covered under the warranty.



Distributed by



APPLETON WOODS LTD

New Lindon House
Catesby Park, Kings Norton
Birmingham B38 8SE
United Kingdom

Sales & Technical Support
Telephone: 0121 458 7740
info@appletonwoods.co.uk

appleton[®]

appleton-scientific.com